QUICK START GUIDE

Moticam 1080 HDMI & USB

EN Quick Start Guide

Thank you for purchasing this Moticam. This camera is a delicate instrument and should be handled with care. When not in use, please store the camera in a safe place with the dust cap attached. Do not let the camera get in contact with water or expose it to extreme temperatures (below 0 or above 50 degrees Celsius). In order to receive the highest level of service, please retain your receipt as proof of purchase and **register your camera online at www.motic.com today.**

Should you have any questions please do not hesitate to **contact our customer service.**

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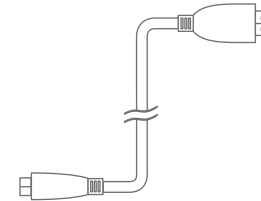
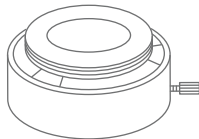
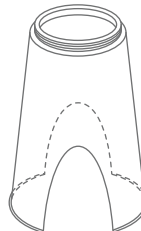
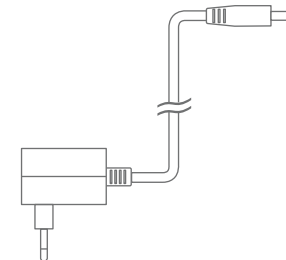
- D**
- EN C-Ring
 - ES Aro-C
 - DE C-Ring
 - FR Anneau type C
 - IT C-Ring

EN Using the camera on an eyepiece

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap from the camera (A) and attach the focusable lens (B).
3. Choose the right size eyepiece adapter (C) and attach this to the focusable lens.
4. Attach this assembly to your microscope's eyepiece by tightening the centering screws on the eyepiece adapter.
5. Connect the USB cable (E) onto the camera and then plug it into the computer.
6. Install the driver if necessary by following the instructions on screen. The driver for the Moticam 1080 is located in the Moticam 1080 Subfolder on the CD.

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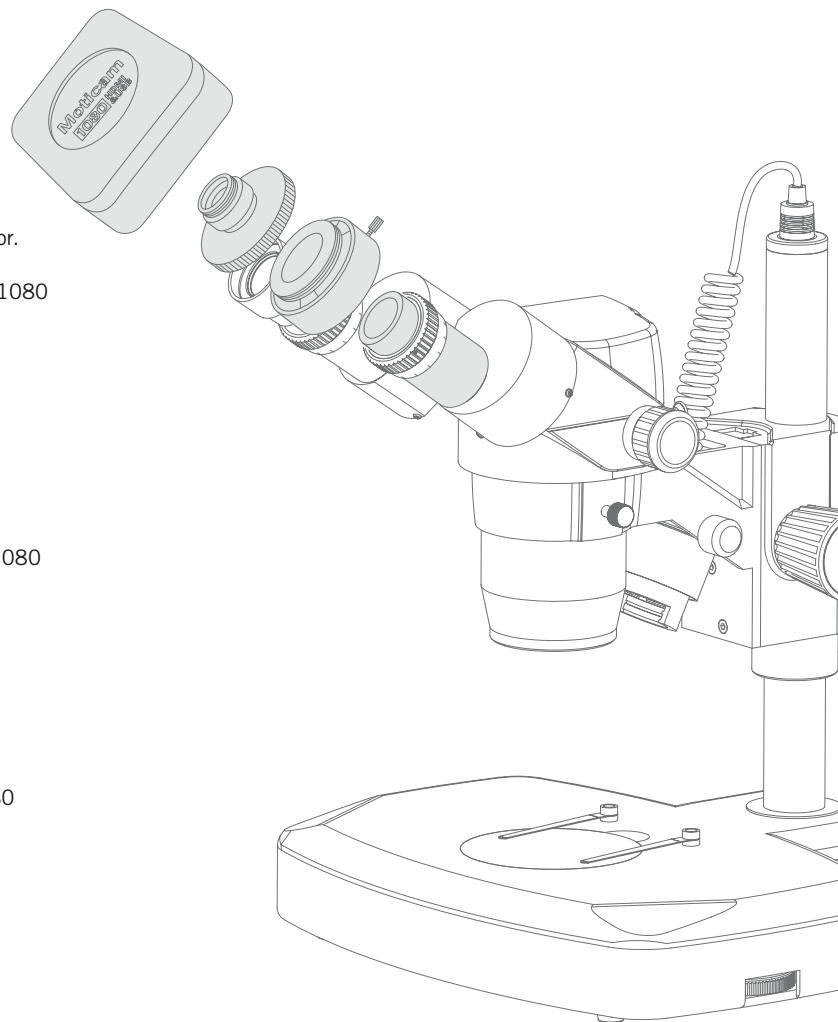
EN Using the camera as a Macro Camera

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap from the camera (A) and attach the focusable lens (B).
3. Attach the Macro Tube (G) onto the lens.
4. Place this assembly over the item you wish to view and ensure there is adequate lighting. If necessary, a desk lamp is a good solution.
5. Connect the USB cable (E) onto the camera and then plug it into the computer.
6. Install the driver if necessary by following the instructions on screen. The driver for the Moticam 1080 is located in the Moticam 1080 Subfolder on the CD.

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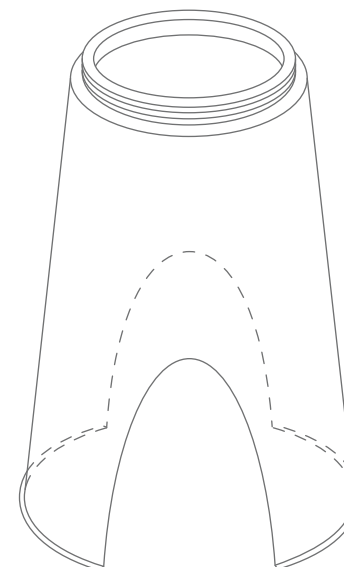
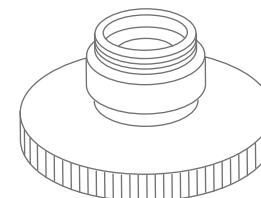
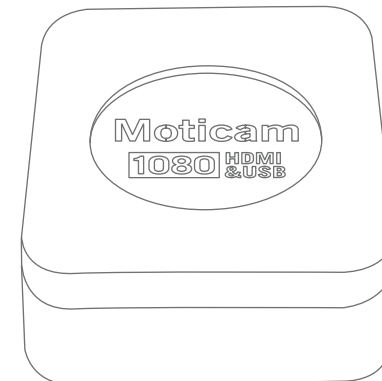
EN Using the camera on a Trinocular Microscope

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap from the camera (A) and attach the microscope's C-Mount adapter (*Please note that the C-Mount adapter is specific to your microscope. For Motic microscopes, please consult the relevant Motic Accessory list and choose an appropriate adapter for your Moticam. Depending on the type of adapter and microscope, the C-Ring (D) may be necessary between the camera and the C-mount adapter).
3. Connect the USB cable (E) onto the camera and then plug it into the computer.
4. Install the driver if necessary by following the instructions on screen. The driver for the Moticam 1080 is located in the Moticam 1080 Subfolder on the CD.

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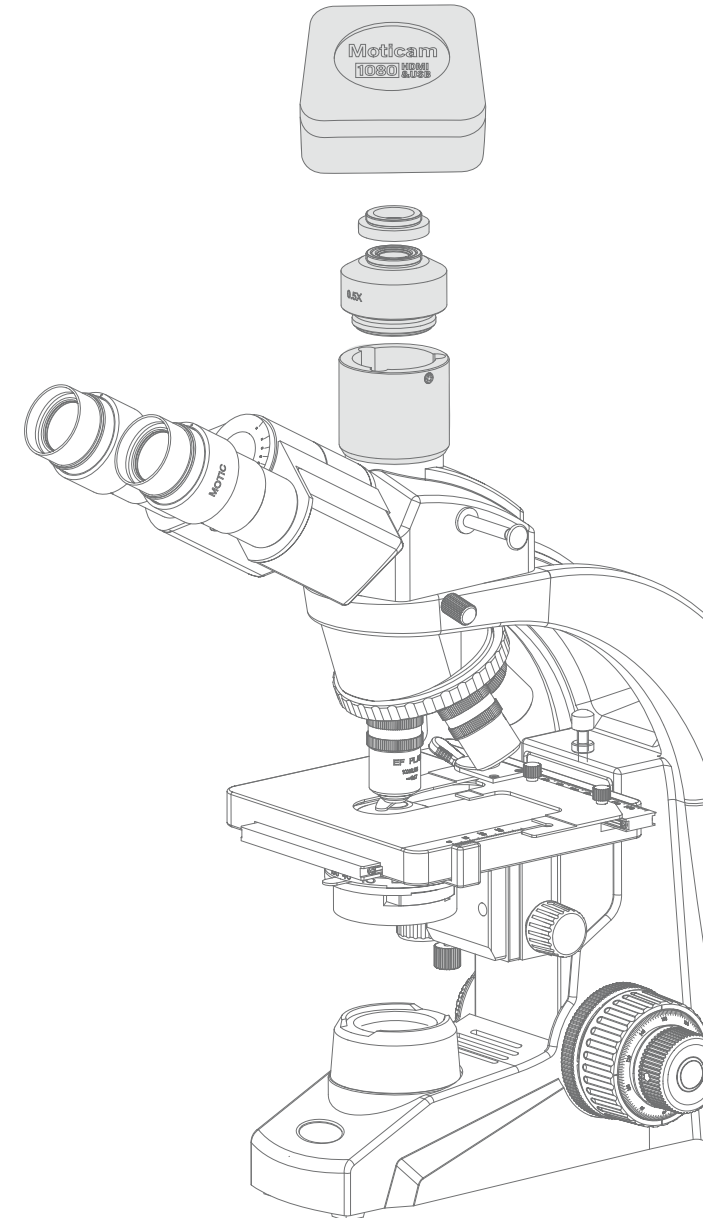
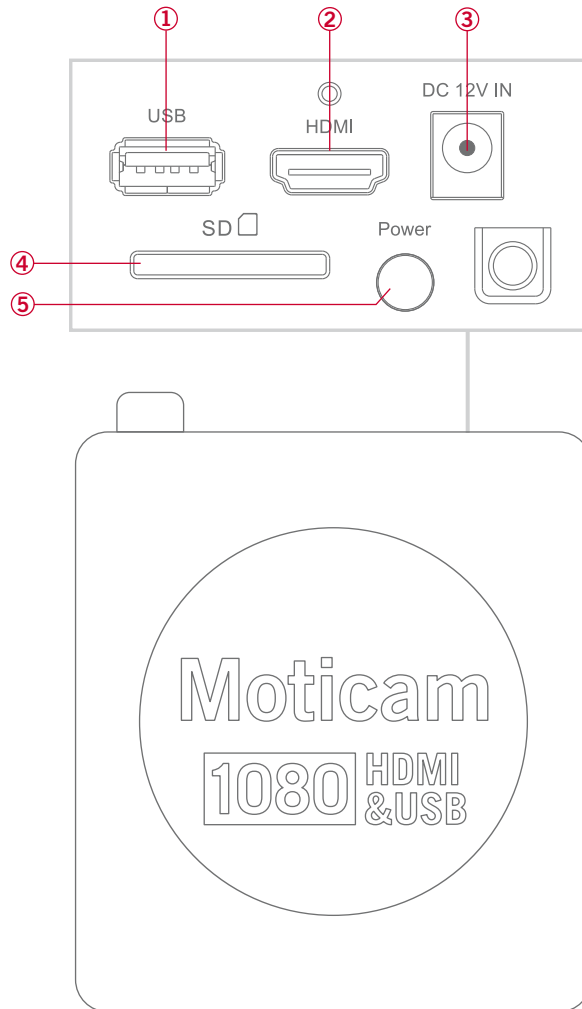
EN 1. The USB connection can be used to connect the Moticam 1080 directly to a computer and use the camera with the provided Motic Images Plus software. This USB port can also be used to connect a USB mouse so that the camera can be controlled without a computer.

2. HDMI port for connecting to an HDMI ready display (the cable is provided in the Moticam 1080 box).

3. Power port for connecting the included Power Supply.

4. SD Card slot for capturing images and video files onto an SD card. The maximum size compatible is 32GB SDHC.

5. Pressing this button will turn the camera on and off.

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EN On-Screen Display

When the camera is connected to an HDMI display and controlled through a USB mouse, you can simply move the mouse cursor to the right to activate the on-screen function display for capturing images and also for adjusting various aspects of the image parameters.

You can also use this setting to change the on-screen display language from the default English to Japanese or Chinese.

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